Quanti-Tray^{*} Sealer

Model 2X • User Manual

For technical support, please call:

North/South America: +1 207 556 4496 or +1 800 321 0207

Europe: +00800 4339 9111

UK: +44 (0) 1638 676800

China: +86 21 61279528

Japan: +81 422 71 5921

Australia: +1300 44 33 99

idexx.com/water

*Quanti-Tray, Defined Substrate Technology, DST, Colilert, Colisure, Enterolert and Pseudalert are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

© 2013 IDEXX Laboratories, Inc. All rights reserved. • 06-03128-04

IDEXX

One IDEXX Drive
Westbrook, Maine 04092 USA

Quanti-Tray* Sealer

Model 2X • User Manual

English

CAUTION: RISK OF ELECTRICAL SHOCK

Do not open the cover of this Sealer.

Do not allow water to spill into this Sealer.

Pour water samples into Quanti-Trays as described in the Quanti-Tray insert.

CAUTION: RISK OF SKIN BURNS

Your company/institution may qualify personnel for cleaning the inside of the Sealer. Qualified personnel should refer to the preventive maintenance instructions for directions on how to clean the inside of the Sealer.

Description

The IDEXX Quanti-Tray* Sealer Model 2X is a motor-driven, heated roller instrument designed to seal IDEXX Quanti-Trays (including Quanti-Tray*/2000). This Sealer, used with Quanti-Trays and any IDEXX reagent in 100mL format, such as Colilert*, Colilert-18*, Colisure*, Enterolert*, Pseudalert*, and HPC for Quanti-Tray*, automates the sample handling of bacterial enumeration. Together, they produce counts as accurate as membrane filtration with the ease of a presence/absence test.

Diagram 1


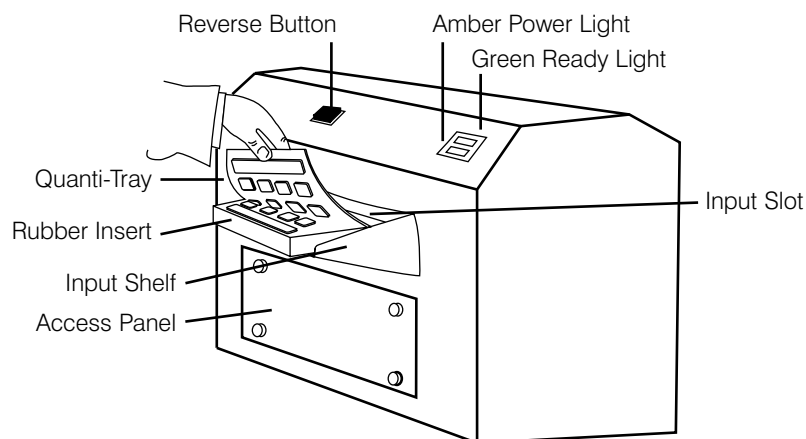
Diagram 2

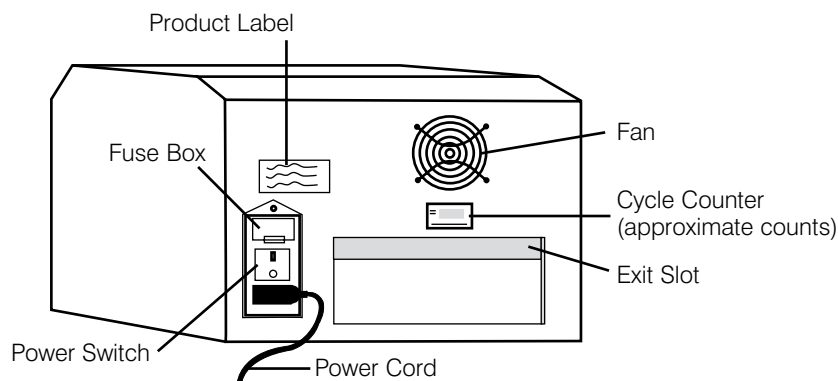Contents

Each Model 2X Sealer comes with the base Sealer unit, an input shelf, one Quanti-Tray* rubber insert, one Quanti-Tray/2000 rubber insert, Most Probable Number (MPN) tables, this user manual and preventive maintenance instructions. The 115-V model also comes with a standard U.S. three-prong grounding power cord. The 230-V model also comes with three cords: a U.K. cord, a European/Shuko cord and a Swiss cord.

Note: Extra Quanti-Tray rubber inserts (catalog number WQTSRBR-51) and Quanti-Tray/2000 rubber inserts (catalog number WQTSRBR-2K) are available separately.

Setting Up

- Unpack contents and save packaging for future shipping.
- Position Sealer on a level surface with adequate space for trays to be inserted and ejected. Sealer can be used in any orientation (left to right or right to left).
- Attach input shelf to Sealer by inserting shelf tabs into the two slots on front of Sealer and clipping firmly in place.
- Ensure that power switch is turned off.
- Plug power cord into Sealer and then into a grounded outlet of proper voltage (see product label).

Operating Directions

WARNING: Personal injury, Sealer damage, property damage and/or inaccurate test results may occur unless Sealer is used according to these instructions.

- Turn power switch on. The amber power light should illuminate.
- Allow the Sealer to warm up and the green ready light to illuminate (up to 20 minutes). Sealer will not operate until both the amber power light and the green ready light are illuminated, indicating that the unit has reached operating temperature.
- Place an empty Quanti-Tray or Quanti-Tray/2000 rubber insert on the input shelf with the large cutout facing away from the Sealer.
- Place a Quanti-Tray or Quanti-Tray/2000 filled with sample and Defined Substrate Technology (DST*) reagent onto the rubber insert, making sure that the tray is properly seated in the rubber insert, with each well of the tray in its corresponding rubber insert hole. **CAUTION: Do not apply labels of any kind to trays prior to sealing. This could cause damage to the Sealer. Use a felt tip marker, or apply label after the tray has been sealed.**
- Slide the rubber insert with tray into the Sealer until the motor grabs the rubber insert and begins to draw it into the Sealer.
- In approximately 15 seconds, the tray will be sealed and partially ejected from the rear of the Sealer. Remove the rubber insert and tray from the rear of the Sealer **immediately**.
- If at any time you wish to reverse the motor drawing the rubber insert into the Sealer (for example, if a misaligned tray is accidentally fed into the Sealer), press and hold the reverse button. However, do not reverse the motor once the rubber insert has been drawn fully into the input slot.
- Multiple rubber inserts can be run consecutively without pausing.
- Turn off Sealer when not in use.

Technical Specifications

| Weight | Dimensions | Ambient Temperature | Power |
|--------|-----------------------------|---------------------|---|
| 35 lb | 12" H x 11" D x 16" W | 32–90°F | 115 V, 60 Hz, 6 amp (part number 99-10893-00) |
| 16 kg | 30 cm H x 27 cm D x 39 cm W | 0–32°C | 230 V, 50 Hz, 3 amp (part number 99-10896-00) |

Cleaning

The rubber inserts may be autoclaved or cleaned with isopropyl alcohol or household bleach, taking the usual precautions when handling such liquids. Be sure to clean the rubber insert if it shows any signs of built-up grime.

Clean the outside of the Sealer with a soft, dry cloth. A soft cloth moistened with water, household bleach, or isopropyl alcohol may also be used, taking the usual precautions when handling such liquids.

The access panel should only be opened by personnel qualified to clean the inside of the Sealer. Refer to the preventive maintenance instructions for directions on how to clean the inside of the Sealer. **Cleaning should only be performed by trained personnel at your facility.** Do not open the access panel or tilt the Sealer if sample is dripping from the unit.

Changing Fuses

Fuses are located in the fuse holder just above the power switch. If it is necessary to change a fuse, turn off the Sealer and unplug the power cord from the Sealer before opening fuse holder. Use 6-amp Bussman MDL-6 fuses or equivalent in the 115-V unit and 4-amp Bussman GDC-4 fuses or equivalent in the 230-V unit.

Troubleshooting

| Symptom | Recommended Action |
|---|--|
| Amber power light doesn't illuminate. | <ul style="list-style-type: none">• Verify that the Sealer is plugged into a live outlet.• Ensure that the power switch is on.• Check fuses and replace if necessary. |
| Amber power light comes on, but the green ready light takes more than 20 minutes to turn on or goes off during operation. | <ul style="list-style-type: none">• Call IDEXX Technical Support or your local distributor. |
| Motor won't start when a Quanti-Tray is inserted. | <ul style="list-style-type: none">• Be sure that Sealer has warmed up and that the green ready light is illuminated.• Be sure that the Quanti-Tray is face-down in the rubber insert with the white tray backing facing upward. |
| Motor starts when Quanti-Tray is inserted but doesn't pull tray through. | <ul style="list-style-type: none">• Be sure tray is in the rubber insert.• Ensure that rubber insert is pushed firmly into roller until motor engages it.• Clean rubber insert. |
| Quanti-Tray and rubber insert get stuck in the Sealer. | <ul style="list-style-type: none">• Press and hold the reverse button to reverse the motor until the tray is ejected from the input slot. Note: If the tray has gone all the way into the Sealer, do not use the reverse button because the tray may become lodged in the Sealer. Instead, pull the rubber insert out of the exit slot. |
| Sealer makes loud hissing noise when sealing, or liquid is dripping from the bottom of the Sealer. | <ul style="list-style-type: none">• This is part of normal operation when trays are overfilled. Do not fill trays with more than 100 mL of sample. The Sealer is designed to automatically remove excess sample from the tray and discard it into the bottom of the Sealer. |
| Trays appear to have blistered, or paper backing is yellowed, after sealing. | <ul style="list-style-type: none">• Call IDEXX Technical Support or your local distributor. Sealer heated roller may be running too hot. |

LIMITED WARRANTY

IDEXX Laboratories, Inc. ("IDEXX") warrants this product to conform to our published specifications, when stored under appropriate conditions and given normal, proper and intended usage, until the expiration of its stated shelf life, or, if none is stated, for one year from the date of delivery of this product to the original end user purchaser ("Buyer"). IDEXX agrees during the applicable warranty period to replace all nonconforming products within 30 days after date of return to IDEXX and without cost to Buyer. IDEXX shall not have any obligation under this Limited Warranty to make replacements which result, in whole or in part, from catastrophe, fault or negligence of the Buyer, or anyone claiming through or on behalf of the Buyer, or from improper use of the products, or use of the products in a manner for which they were not designed, or by causes external to the products.

Buyer shall notify IDEXX of any products which it believes to be nonconforming during the warranty period. At IDEXX's option, such products shall be returned by Buyer, transportation and insurance prepaid, to IDEXX's designated facility for examination and testing. IDEXX shall repair or replace, within 30 days of receipt by IDEXX, any such product found to be so nonconforming and promptly return such products to Buyer, transportation and insurance prepaid. Should IDEXX's examination and testing not disclose any nonconformity covered by the foregoing warranty, IDEXX shall so advise Buyer and dispose of or return the product in accordance with Buyer's instructions and at Buyer's sole expense.

THE PROVISIONS OF THE FOREGOING LIMITED WARRANTY ARE IN LIEU OF ANY OTHER WARRANTY, WHETHER EXPRESS OR IMPLIED, WRITTEN OR ORAL (INCLUDING ANY WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE). IDEXX'S LIABILITY ARISING OUT OF THE MANUFACTURE, SALE OR SUPPLYING OF THE PRODUCTS OR THEIR USE OR DISPOSITION, WHETHER BASED UPON WARRANTY, CONTRACT, TORT OR OTHERWISE, SHALL NOT EXCEED THE ACTUAL PURCHASE PRICE PAID BY BUYER FOR THE PRODUCTS. IN NO EVENT SHALL IDEXX BE LIABLE TO BUYER OR ANY OTHER PERSON OR ENTITY FOR SPECIAL, INCIDENTAL, CONSEQUENTIAL, INDIRECT OR EXEMPLARY DAMAGES (INCLUDING, BUT NOT LIMITED TO, LOSS OF PROFITS OR LOSS OF USE DAMAGES) ARISING OUT OF THE MANUFACTURE, SALE OR SUPPLY OF THE PRODUCTS. THE FOREGOING WARRANTIES EXTEND TO BUYER ONLY AND SHALL NOT BE APPLICABLE TO ANY OTHER PERSON OR ENTITY INCLUDING, WITHOUT LIMITATION, CUSTOMERS OF BUYER.

For technical support, please call:

North/South America: +1 207 556 4496 or +1 800 321 0207

Europe: +00800 4339 9111

UK: +44 (0) 1638 676800

China: +86 21 61279528

Japan: +81 422 71 5921

Australia: +1300 44 33 99

idexx.com/water

*Quanti-Tray, Defined Substrate Technology, DST, Colliert, Colisure, Enterolert and Pseudalert are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

© 2013 IDEXX Laboratories, Inc. All rights reserved. • 06-03128-04

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© 2014 IDEXX Laboratories, Inc. Todos os direitos reservados. • 06-03128-04