

## IDEXX xChek\*: Creating a User-Defined Assay for the IDEXX Influenza Virus Antibody Test - US and Canada

Creating a user-defined assay is an easy process of two tasks—editing an xChek\* file to allow you to create an assay, and then creating the assay within xChek software.


### To edit the xChek.ini file:

1. From the task bar, choose **Start> Run**, and type “xchek.ini” into the text box. Click **OK**.
2. After the xChek.ini file opens, locate the line labeled “AllowNewAssays.”
3. If the value for that line is “0,” change it to “1.” If the value is already “1,” do not do anything.
4. Locate the section labeled “[BasicGraphs],” and scroll to the end of the section.
5. Enter “SIV=17.”
6. From the menu bar, choose **File> Save**, and then choose **File> Exit**.

### To create the assay:


1. From the task bar, choose **Start> Programs> xChek> xChek** to open the xChek program.
2. If prompted to log on, enter your initials and then click **OK**.
3. From the menu bar, choose **Database> Assays> All** to open the Assay dialog box.
4. Click the **New** button, and enter “Influenza Virus Antibody Test - US and Canada” into the **Name** field, “SIV” into the **Code** field, and “unknown” into the **Kit Lot** field. Today’s date (mm/dd/yyyy) is automatically entered into the Expiration field.

**NOTE:** When you begin testing, update the information in the Kit Lot and Expiration fields with the actual kit lot number and expiration date.

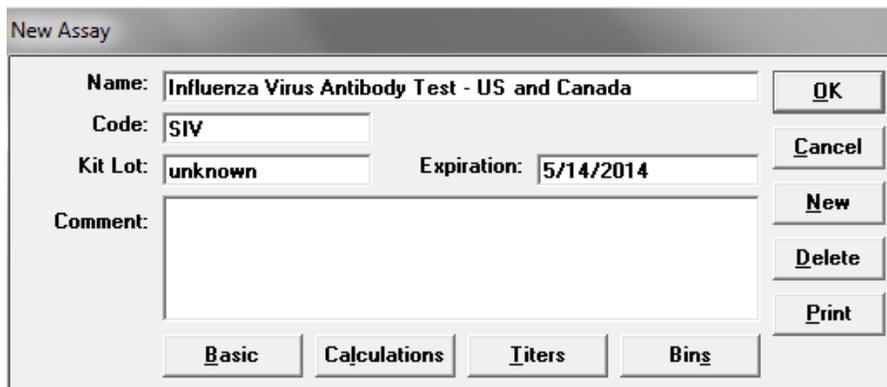

5. Click the **Basic** button, and select or enter the following information in the respective fields:

Case Type: Animal	Samp Filter: 650
Template: SIV Vert	Ref. Filter: 0
Species: Cattle	Dilution: 10
Technology: ELISA	Wells: 1

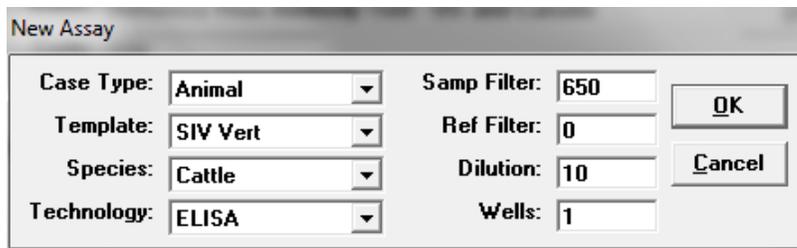

**Note:** Because of a problem with selecting swine as the species type in versions 3.2 and 3.3 of xChek software, the species used for this user-defined assay is cattle.

- Click **OK** to save the settings.
- Click the **Calculations** button, and select or enter the following information:

Formulas:

"S" Part of Ratio: Sample1

"P" Part of Ratio: None

"N" Part of Ratio: Negative

Blocking Factor: None

First Calculation:

Variable: S/N

Positive Cutoff: 0.6

Suspect Cutoff: 0.6

Comparison: <

Second Calculation:

Variable: None

Positive Cutoff: 0

Suspect Cutoff: 0

Comparison: None

**New Assay**

**Formulas**

"S" Part of Ratio: Sample1

"P" Part of Ratio: None

"N" Part of Ratio: Negative

Blocking Factor: None

**First Calculation**

Variable: S/N

Positive Cutoff: 0.6

Suspect Cutoff: 0.6

Comparison: <

**Second Calculation**

Variable: None

Positive Cutoff: 0

Suspect Cutoff: 0

Comparison: None

OK  
Cancel  
Reset


- Click **OK** to save these settings, and then click **OK** to save the assay to the database. You can now test for influenza antibodies using the xChek Assay Management System.

**NOTE:** xChek does not evaluate controls or results for validity when you use a user-defined assay. You must evaluate the results from each assay in accordance with good laboratory practices. To evaluate your assay validity, refer to the insert provided with the test.


### Setting Up Additional Display Variables

Only OD values are displayed in the reporting options for this user-defined assay. To set up additional display variables, follow the directions below.

- From the xChek menu bar, choose **Reports> Analyze Cases** to open the Filter Criteria for Analyze dialog box.
- Enter the desired search criteria and click **OK**. The Select Cases for Analyze dialog box appears.
- Select the desired cases and click **OK**. The Analyze Report dialog box appears.

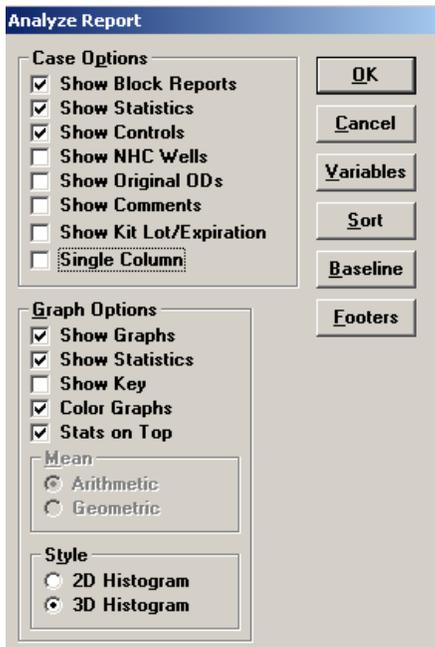

4. Click the **Variables** button to open the Display Variables dialog box.
5. Select the desired options and click **OK**. Click **OK** again to save the settings.

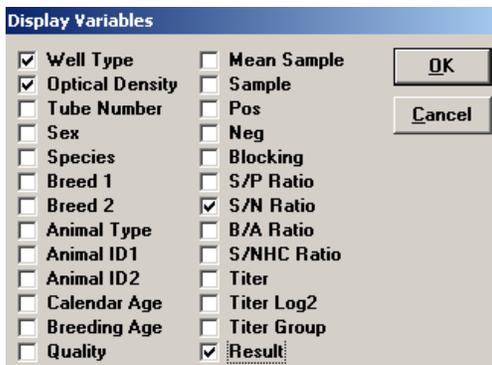

For more information, contact IDEXX Technical Services at 1-800-548-9997, option 2, or visit us on the web at [idexx.com](http://idexx.com).

Test With Confidence™

**IDEXX**

One IDEXX Drive  
Westbrook, Maine 04092 USA

© 2014 IDEXX Laboratories, Inc. All rights reserved. • 105262-00

\*xChek and Test With Confidence are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries. The IDEXX privacy policy is available at [idexx.com](http://idexx.com).